

# Heart Weight RBRCLA\_HWT\_001

## Purpose

To evaluate cardiac size using heart weight and body weight.

## Experimental Design

- **Minimum number of animals** : 7M + 7F
- **Age at test:** Week 57

## Equipment

- fine forceps
- surgical scissors
- fine surgical scissor
- kim wipes (tissues) or surgical compress
- laboratory balance
- labelled jar with fixative
- corkplate or wax board
- pins
- jar containing tap water to rinse the tools

## Procedure


Methods and procedures used not including center-specific data entry methods.

1. Sacrifice the mouse
2. Weigh and record total body weight
3. Place mouse on its back and pin the mouse onto board with extended extremities (inner side of hands and foot)
4. Wipe or wet the mouse with 70% ethanol to control hair and dander
5. Option A (for mice that are undergoing complete necropsy):
  - Proceed with a complete necropsy and tissue collection according to Centre-specific technical SOP including removal of the heart by dissecting the aortic root immediately above the aortic valves and the superior vena cava above the atria
  - Remove adjacent mediastinal fat pads from the excised heart carefully with forceps
  - Empty heart blood by tapping the heart on a kim wipe (absorbent pad) or surgical compress. Repeat until the heart is totally dry

- Weigh the heart, record the weight in the Centre-specific database, and place the heart in fixative

6. Option B (for mice that are not undergoing complete necropsy):

- Open the skin in the ventral midline and in the direction of the extremities and extend cut to hands and feet
- Open the muscular abdominal wall in the midline and along the lower margin of the rib cage with small forceps
- Open rib cage by removing the sternum and adjacent ribs
- Remove the heart by dissecting the aortic root immediately above the aortic valves and the superior vena cava above the atria
- Remove adjacent mediastinal fat pads from the excised heart carefully with forceps
- Empty heart blood by tapping the heart on a kim wipe (absorbent pad) or surgical compress. Repeat until the heart is totally dry
- Weigh the heart and record the weight in the Centre-specific database
- Discard the heart

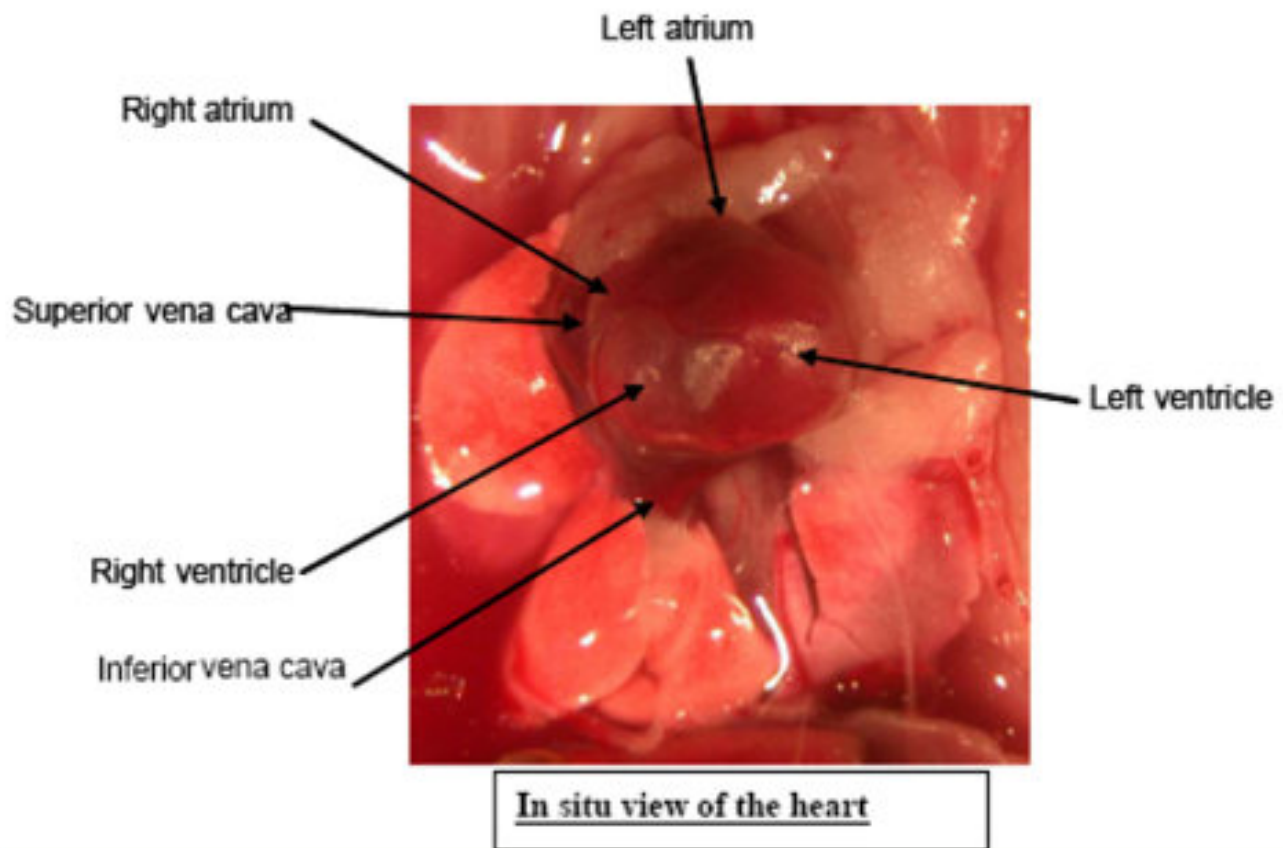

## Notes

All data are collected at a local workstation in the necropsy room (attached to a digital balance) and uploaded to the Centre-specific pathology data capture system.

### Data QC

Mouse weight between 5 grams and 150 grams

## IMPC Parameters (+ontology annotations)

Weight (in mg)

## Data Analysis, annotation and display (+statistics)

Statistics: ANOVA/Wilcoxon test using normalized heart weights (for body weight) displayed in boxplot

## Parameters and Metadata

### Heart weight RBRCLA\_HWT\_008\_001 | v1.2

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Unit Measured: mg

---

### Body weight RBRCLA\_HWT\_007\_001 | v1.3

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: false

Unit Measured: g

---

### Heart weight normalised against tibia length RBRCLA\_HWT\_013\_001 | v1.3

simpleParameter

Req. Analysis: false

Req. Upload: false

Is Annotated: false

Derivation:  $\text{div}(\text{'RBRCLA\_HWT\_008\_001'}, \text{'RBRCLA\_HWT\_002\_001'})$

Date equipment last calibrated RBRCLA\_HWT\_009\_001 | v1.2

procedureMetadata

Req. Analysis: false

Req. Upload: false

Is Annotated: false

Equipment ID RBRCLA\_HWT\_006\_001 | v1.0

procedureMetadata

Req. Analysis: false

Req. Upload: true

Is Annotated: false

Heart weight normalised against body weight RBRCLA\_HWT\_012\_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: false

Is Annotated: false

Derivation:  $\text{div}(\text{'RBRCLA\_HWT\_008\_001'}, \text{'RBRCLA\_HWT\_007\_001'})$

# Equipment model RBRCLA\_HWT\_011\_001 | v1.0

procedureMetadata

Req. Analysis: true                      Req. Upload: true                      Is Annotated: false

Options: PG3001-S, Scout Pro SPU123, MS104S, QUINTIX124-1S, Adventurer AX223/E, TP-114, P-403, AV213C, AV212C, Adventurer Pro, AV2101, AB104-S, GF-200, HR-120, TE212, 201-10, BCE124I-1SJP,

-----

# Experimenter ID RBRCLA\_HWT\_003\_001 | v1.0

procedureMetadata

Req. Analysis: false                      Req. Upload: true                      Is Annotated: false

# Equipment manufacturer RBRCLA\_HWT\_010\_001 | v1.0

procedureMetadata

Req. Analysis: true                      Req. Upload: true                      Is Annotated: false

Options: Denver Instrument, A & D, Mettler Toledo, Sartorius AG Germany, Ohaus, Radwag,

-----

# Tibia length RBRCLA\_HWT\_002\_001 | v1.1

simpleParameter

Req. Analysis: false                      Req. Upload: false                      Is Annotated: true

Unit Measured: mm

---

**Method of sacrifice** RBRCLA\_HWT\_005\_001 | v1.1

procedureMetadata

Req. Analysis: false      Req. Upload: true      Is Annotated: false

**Options:** Ketamine (137mg/kg)/Xylazine (6.6mg/kg),  
Ketamine (100mg/kg)/Xylazine (10mg/kg), Carbon dioxide, Isoflurane overdose,  
Ketamine (110mg/kg)/Xylazine (11mg/kg), None, Avertin, Cervical dislocation,  
Cardiac puncture, Anesthetized, Pentobarb (0.1ml),  
Ketamine (110mg/kg)/Xylazine (11mg/kg)/ Antisedan (1mg/kg),  
Ketamine(100mg/kg)/ Xylazine (10mg/kg )/ Antisedan (1mg/kg), Exsanguination,  
Medetomidine/Midazolam/Butorphanol,

---

**Date of sacrifice** RBRCLA\_HWT\_001\_001 | v1.0

procedureMetadata

Req. Analysis: false      Req. Upload: true      Is Annotated: false

---